



Induced Pluripotent Stem Cells (Control Lines from CD34+ Blood Cells; Female)

Product Information

Catalog Number	ASE-9110 (Female)
Description	Applied StemCell, Inc. provides fully characterized control iPSC lines that are derived using a standardized integration-free, episomal reprogramming protocols from CD34+ cord blood cells. The ASE-9110 control line is from a female donor, and has been characterized for the expression of pluripotency markers. They are ideal as isogenic controls for disease modeling, drug and toxicity screening.
Tissue	CD34+ Cord Blood Cells
Sex	Female
Race	Caucasian
Clinical Information	Normal
Quantity	≥1.0x10 ⁶ viable cells/vial
Shipping	Dry ice
Storage and Stability	Store in liquid nitrogen freezer immediately upon receipt. This product is stable for at least 6 months from the date of receiving when stored as directed.
Quality Control	Each lot of human iPS cells has been tested for growth and viability following recovery from cryopreservation. In addition, the iPSCs have also been tested for normal female karyotype, expression of stem cell markers (SOX2, NANOG, OCT4, TRA-1-81, and TRA-1-60; Figure 1) and for the absence of mycoplasma and pathogens (CoA available upon request).
Safety Precaution	PLEASE READ BEFORE HANDLING ANY FROZEN VIALS. Please wear the appropriate Personal Protection Equipment (lab coat, thermal gloves, safety goggles and a face shield) when handling the cells. Handle the frozen vials with due caution. Please be aware that the following scenario can occur: Liquid nitrogen can leak into the vials when the vials are submerged in liquid nitrogen. Upon thawing, the liquid nitrogen returns to the gas phase, resulting in a dangerous build-up of pressure within the vial. This can result in the vial exploding and expelling not only the vial contents but also the vial cap and plastic fragments of the vial.
Restricted Use	This product is for research use only and not intended for human or animal diagnostic or therapeutic uses.

Applied StemCell, Inc.

521 Cottonwood Dr. #111, Milpitas, CA 95035

Phone: 866-497-4180 (US Toll Free); 408-773-8007 Fax: 408-773-8238

info@appliedstemcell.com www.appliedstemcell.com

Copyright 2019, Applied StemCell, Inc. All rights reserved. This information is subject to change without notice.

Characterization of Control iPSC Lines

Expression of Pluripotency Markers

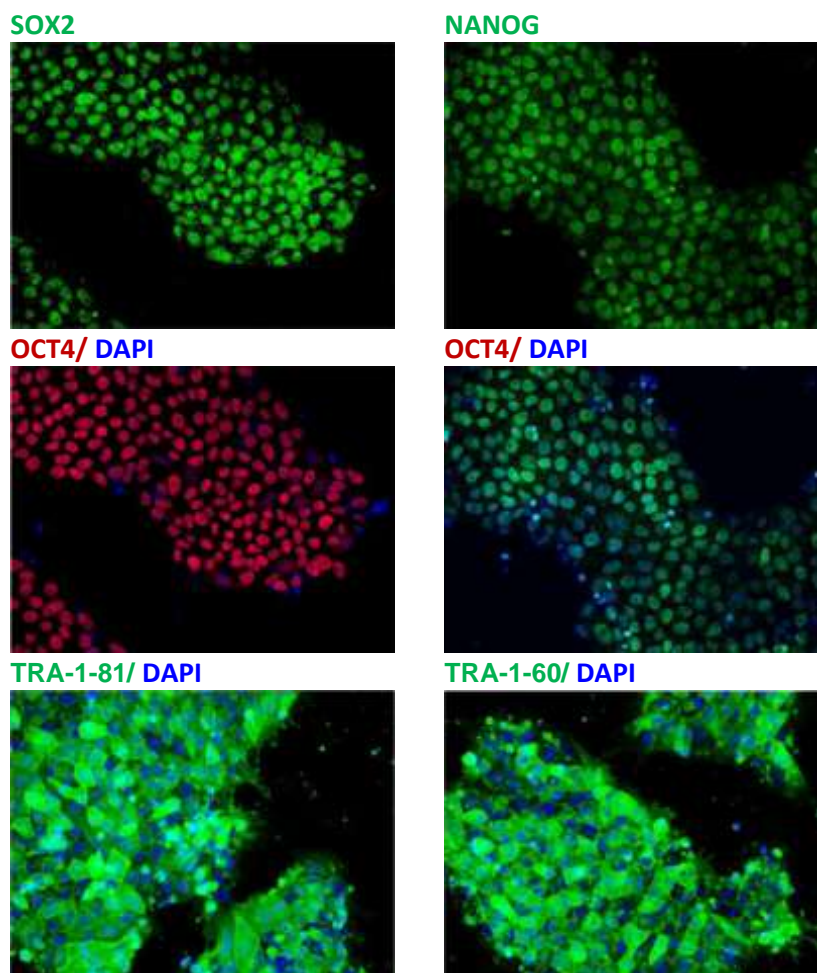


Figure 1. Expression of pluripotency markers in Control iPSC line, ASE-9110. The control iPSC line, derived from female cord blood cells expresses pluripotency markers NANOG, OCT4, SOX2, TRA-1-60, and TRA-1-81, indicating pluripotency of the iPSC line. Nucleus stained with DAPI (blue).

Additional Reagents Required but not Provided

- Rock Inhibitor, Stemgent, Cat# 04-0012
- mTeSR™1 Feeder Free Media, STEMCELL Technologies, Cat# 85850
- Corning® Matrigel® hESC-Qualified Matrix, *LDEV-free, Corning, Cat# 354277
- DMEM, high glucose, pyruvate, ThermoFisher, Cat# 11995081
- StemPro™ Accutase™ Cell Dissociation Reagent, ThermoFisher, Cat# A1110501
- CryoStem™ Freezing Medium, Stemgent, Cat# 01-0013-50

Equipment Required

Equipment	Specifications
Centrifuge	(e.g. Thermo IEC Centra CL2)

Conical Tubes	Polystyrene Conical Tube 15 mL, Corning Inc., Cat # 430790 Polystyrene Conical Tube 50 mL, Corning Inc., Cat# 176740
Cell Counter/Hemocytometer	
Cell Culture Dishes	60mm/15mm, Corning Inc., Cat# 430196

Protocol

All steps must be performed according to standards required for sterile cell culture work.

1. Coating Cell Culture Dishes or Flasks with Corning® Matrigel® hESC-qualified Matrix Substrate

Please read the producer' manual for handling of Corning® Matrigel® hESC-qualified Matrix.

Important producer's notes:

It is extremely important that Corning® Matrigel® hESC- qualified Matrix and all culture ware or media coming in contact with Corning® Matrigel® hESC-qualified Matrix should be pre-chilled/ice-cold since Corning® Matrigel® hESC-qualified Matrix will start to gel above 10°C.

The dilution is calculated for each lot based on the protein concentration. Prepare aliquots according to the dilution factor provided on the Certificate of Analysis (use 1.5-2.0 mL tubes). The volume of the aliquots is typically between 270-350 µL.

- 1.1 Pre-chill pipette tips and dishes at 4°C.
- 1.2 Thaw an aliquot (typically between 270-350 µL) of Corning® Matrigel® hESC-qualified Matrix at 4°C (approximately 45-60 minutes).
- 1.3 Transfer the aliquot on ice into a biological safety cabinet.
- 1.4 Prepare a 25 mL aliquot of cold DMEM in 50 mL conical tube and keep on ice.
- 1.5 Using a p1000 micropipette, transfer 1000 µL of the cold DMEM from the above tube into the tube with Matrigel, and mix up and down several times. Transfer the Matrigel solution to the 50 mL conical tube containing cold DMEM and mix several times with a serological pipette (keep on ice).
- 1.6 Immediately, coat pre-chilled culture dishes with Matrigel/DMEM solution (volume 120-150 µL/cm²).
- 1.7 Distribute coating matrix evenly and incubate at room temperature (15-25°C) for at least 1 hour before use.
- 1.8 Coated dishes can be used immediately or can be stored at 4°C for up to 7 days (aseptic conditions).

2. Recovering iPSC from Frozen Stock (Feeder-free, Single Cell Suspension)

- 3.1 Pre-warm mTeSR1 culture medium to 37°C, and supplement the medium with 10 µM (final concentration of ROCK inhibitor).
- 3.2 Pre-warm Matrigel coated dishes.
- 3.3 Pipette out 5 mL of the above medium into a 15 mL conical tube.
- 3.4 Transfer a vial with frozen iPSC (on dry ice) to the operation side and thaw cells immediately in 37°C degree water bath, until only a small piece of ice is still visible (approximately 1-1.5 minutes).
- 3.5 Transfer the vial to the biological safety cabinet and transfer thawed iPSC slowly and drop-wise into a 15 mL conical tube containing 5 mL of pre-warmed mTeSR1 + ROCK medium under constant swirling.
- 3.6 Wash the cryo-vial with additional 1 mL of medium and transfer to the same 15 mL tube.
- 3.7 Centrifuge cells at 1000 rpm for 5 minutes.
- 3.8 Aspirate supernatant carefully without disturbing the pelleted cells and re-suspend the pellet carefully but thoroughly (5-6 up-and-down mixes) in 5 mL of fresh mTeSR1 + ROCK medium using 5 mL serological or p1000 micro-pipette with a long tip.
- 3.9 Count the cells (optional).
- 3.10 Aspirate Matrigel solution from coated dishes and immediately seed iPSC at 80,000-150,000 cells/cm² (do not allow the matrix to dry out). We recommend seeding the entire vial onto one 60 mm dish.

Optional: Aspirate Matrigel before the cells are thawed and add enough medium to cover the bottom of the culture dish.

- 3.11 Distribute cells evenly by performing cross-like moves and place dish(es) in 37°C / 5%CO₂ cell culture incubator.
- 3.12 Change mTeSR1 medium (without ROCK inhibitor) every day.

3. Enzymatic Passage of Feeder-Free Cultured iPSC Using Accutase (Single Cell Suspension)

- 4.1 Pre-warm plates coated with Matrigel at 37°C.
- 4.2 Prepare new plate for transfer of clones: Remove the Matrigel solution from the plates and add appropriate volume of pre-warmed mTeSR1 supplemented with 10 µM ROCK inhibitor to each plate (5 mL for 60mm, 10 mL for 100mm plates, etc.).
- 4.3 Aspirate media from the plate to be passaged. Add pre-warmed Accutase to each plate (2.5 mL for 60mm, 5 mL for 100mm plates), and place the plate in a 37°C incubator. Observe intermittently to determine when cells begin to detach. Do not leave cells in Accutase for longer than 10 minutes.
- 4.4 When cells are detached, add an equal volume of pre-warmed DMEM to each plate to dilute the Accutase (2.5 mL for 60mm, 5 mL for 100mm plates).
- 4.5 Pipette the cell mixture in the plates up and down using a p1000 pipette. Wash the bottom of the plate well to ensure detachment of cells. Transfer the cell suspension to a conical tube.
- 4.6 Centrifuge tubes at 1000 rpm for 5 minutes.
- 4.7 Aspirate the supernatant and re-suspend cells thoroughly using p1000 pipette in mTeSR1 (4 mL for pellet from 60mm dish, 9 mL for pellet from 100mm dish).
- 4.8 Centrifuge tube at 1000 rpm for 5 minutes.
- 4.9 Take 10 µL cell mixture, mix with 10 µL of Trypan blue and use for cell count.
- 4.10 Plate the cells on the previously prepared Matrigel coated plates with mTeSR1 + ROCK medium at densities ranging from 50,000-100,000 cells/cm².
- 4.11 Incubate at 37°C/ 5%CO₂, and observe attachment within 12 hours. Change media after 24 hours and daily, thereafter.

4. Cryopreservation of iPSCs

- 4.1 Aspirate media from plates to be harvested. Add pre-warmed Accutase to each plate and place the plate in a 37°C incubator. Observe intermittently to determine when cells begin to detach (approximately 5 minutes). Do not leave cells in Accutase for longer than 10 minutes.
- 4.2 When cells are detached, add an equal volume of pre-warmed DMEM to each plate to dilute the Accutase, perform cell count and calculate how many vials need to be frozen down (e.g. 2.5 x 10⁶ cells per vial).
- 4.3 Centrifuge the tube at 1000 rpm for 5 minutes.
- 4.4 Remove supernatant from tube and re-suspend cells in appropriate volume of CryoStem freezing media to achieve 2.5 x 10⁶ cells per 1 mL.
- 4.5 Transfer 0.5 mL of the cell suspension in Cryostem™ freezing media to each cryo-vial and place the vials in a freezing container and immediately place the container at -80°C.
- 4.6 Transfer vials to LN2 within a week's time window.